


Electrode to Connector Map	
Assembly type	Deep brain mapping array for acute experiment.
Version	V1.0 Jan 2009.
Connector	Omnetics nano NPD series (A7695-001)
Mating Connector	Omnetics nano NPD series (A7519-001)
Compatible Headstages	TDT RA16CH, LP16CH
	Multi-Channel Systems Two MPA8I
	Plexon HST/16o25-18P, HST/16V-G20
	Neuralynx HS-16
	Alpha Omega 16 mini, TeleSpike
	DataWave Technologies DW-T-HS16, DW-W-HS16, DW-M-HS16
Compatible Adaptors	TDT AC-CH
Note	1 st Site is 0.5mm from the tip. Alcohol sterilization may be incompatible with some of the manufacturing components. Ethylene oxide gas sterilization is recommended.


CONNECTOR PIN OUT

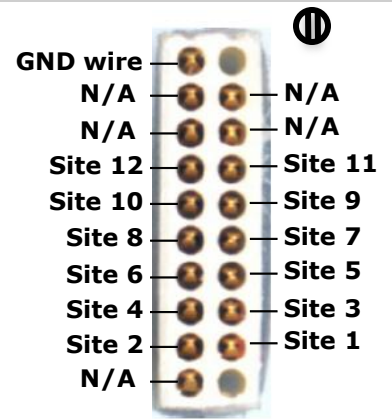

ARRAY PHOTO



ARRAY SITE LAYOUT

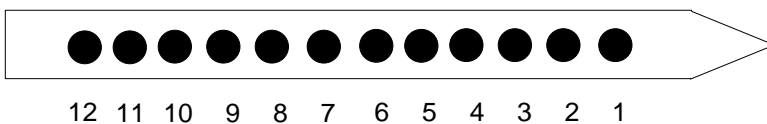

Site-to-site spacing is either 150um or 300um. Please check your order for detail.

